

PRODUCT INFORMATION

MycoDecon

Cat. No.: 34206

PRODUCT DESCRIPTION

Highly effective disinfectant, particularly active against mycoplasmas, but also against bacteria, virus and fungi.

Application: MycoDecon is special spray formulation for the use on any laboratory surfaces and equipment, including benches, incubators, storage boxes of cells, containers for liquid nitrogen.

Decontamination of surfaces:

- Spray MycoDecon directly on the contaminated surface until it is moistened completely.
- After 5 min dry the surface with a clean paper towel.
- **IMPORTANT:** Remove completely, do not leave any residue.

**Please note: May cause discoloration of sensitive surfaces.
Test on small part of the surface prior to use.**

Decontamination of laboratory devices:

- Apply MycoDecon on a paper towel and clean all surfaces of the devices thoroughly.
 - Dry with a clean paper towel.
-

Features

- Ready-to-use
 - Alcohol based
 - Non-corrosive
 - Non-carcinogenic
-

Storage Store at + 15 °C to + 30 °C.

Ver. 07/15